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Revision changes

1. Goal

To describe how feces must be prepared and stored within in the Radboud Biobank, so that:

1. Potential users know how the biomaterial has been handled.
2. The laboratory preparing the biomaterial for storage knows the minimal requirements for the handling of it.

1.1 Scope of application

This procedure is applicable to all employees concerned at the Radboud Biobank and the Radboud Laboratory for Diagnostics.

2. Protocol feces Standard

Collection

Delivery In Feces container.

Type of container Sterile 60 ml PP container, Corning TP35C-013 or equivalent.

Temperature Keep containers at 4°C until preparation.

Preparation

Time until freezing Maximum 72 hours after "collection" if kept refrigerated (4°C).

Temperature 4°C (in refrigerator (4°C) or on ice).

Preparation protocol Homogenize feces by using a wooden spatula. If the feces is too hard to homogenize do not homogenize but divide into portions and note the deviation in the Biobank management system.

Storage

Register Register tubes with feces in the Biobank management system and attach a cryovial label.

Aliquotation 6 aliquots of 1-2 grams should be stored. If there is not enough material store, less aliquots. If the feces is too hard divide the material over the tubes (max. 6, volume may deviate).

Type of tube 7ml polyethylene feces tube, 47x19PC+CAP (Sarstedt 71.9923.610).

Temperature -80°C

Miscellaneous

Deviations Note all deviations from this procedure in the Biobank management system.

3. Quality assurance

Review takes place via an internal audit (see the audit plan).

3.1 Performance indicator

95% of the relevant samples have been prepared and stored within the parameters of this procedure.

4. Comments

NB: This procedure has been written for research focusing DNA analysis of the feces microbiome. If the intended purpose is Metabolomics studies it is imperative that the samples are prepared and frozen as fast as possible with a maximum of 48 hours after collection.

5. Relevant documents

Qdocs, literature, legislation and codes of conduct, website etc.

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